TU-20HT
Thermoregulator
Operator's Manual

Issue 7

Date of issue:

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
3rd May 2005

Techne is a trade mark

© Techne 2002

CONTENTS

TU-20HT Operator's Manual	
	<i>page</i>
SAFETY and INSTALATION	
English	4
Français	6
Deutsch	8
Español	10
THE TU-20C	12
Constant Temperature Baths	12
Description	12
Unpacking	13
Set Up	13
SPECIFICATION	
TU-20HT	18
Baths	19
OPERATION	20
When you switch on the TU-20HT	21
Front Panel controls	21
Setting the Operating Temperature	24
After Use	24
Setting the overtemperature cutout	25
RS232 Serial Interface	26
Sending Data to a PC	26
ADDITIONAL INFORMATION	27
Operator maintenance	27
Calibration	28
Replacement parts	29
Accessories	29

NOTES

Introduction

Please read all the information in this booklet before using the unit.

Warning

HIGH TEMPERATURES ARE DANGEROUS: they can cause serious burns to operators and ignite combustible material.

Techne have taken great care in the design of these units to protect operators from hazards, but Operators should pay attention to the following points:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS;
- DO NOT put hot objects on or near combustible objects;
- DO NOT operate the unit close to inflammable liquids or gases;
- DO NOT place any liquid directly in your unit;
- At all times USE COMMON SENSE.

Operator Safety

All Operators of Techne equipment must have available the relevant literature needed to ensure their safety.

It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures. If the equipment is used in a manner not specified by Techne the protection provided by the equipment to the Operator may be impaired.

All Techne units have been designed to conform to international safety requirements and are fitted with an overtemperature cutout. On some models, the cutout is adjustable and should be set to suit the application. On all other models the cutout is preset to protect the unit.

If a safety problem should be encountered, switch off at the mains socket and remove the plug from the supply.

Installation

1. All Techne units are supplied with a power cable. This may be integral or plug-in.
2. Before connecting the mains supply, check the voltage against the rating plate. The rating plate is on the rear of the unit. Connect the mains cable to a suitable plug according to the table below.

Note that the unit must be earthed to ensure proper electrical safety.

Connections	220V-240V	110V-120V
Live	Brown	Black
Neutral	Blue	White
Earth	Green/yellow	Green

The fused plug supplied with the mains lead for use in the UK is fitted with the following value fuse to protect the cable: 10AMP.

The fuse in the unit protects the unit and the operator

Note that units marked 230V on the rating plate work at 220V; units marked 120V work at 110V. In both cases, however, the heating rate will degrade by approximately 8%.

3. Plug the mains cable into the socket on the rear of the unit.
4. Place the unit on a suitable bench or flat workspace, or in a fume cupboard if required, ensuring that the air inlet vents are free from obstruction.
5. Note that the following symbols may be next to the indicator lamps on the front panel of the units and have the following meanings:

~	: the power indicator
⏏	: the heater indicator
⚡	: the overtemperature indicator

6. Symbols on or near the power switch of the unit have the following meanings:

I	: mains switch On
O	: mains switch Off

After use

When you have finished heating samples, remember that parts of the unit – the tubes, blocks and associated accessories – may be very hot. Take the precautions listed earlier.

Guarantee

The unit is guaranteed against any defect in material or workmanship for the period specified on the enclosed guarantee card. This period is from the date of purchase, and within this period all defective parts will be replaced free of charge provided that the defect is not the result of misuse, accident or negligence. Servicing under this guarantee should be obtained from the supplier.

Notwithstanding the description and specification(s) of the units contained in the Operator's Manual, Techne hereby reserves the right to make such changes as it sees fit to the units or to any component of the units.

This Manual has been prepared solely for the convenience of Techne customers and nothing in this Instruction Book shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the units or components.

Operator maintenance

NOTE: THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL.

REMOVING THE SIDE, FRONT OR REAR PANELS EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES.

THERE ARE NO OPERATOR MAINTAINABLE PARTS WITHIN THE EQUIPMENT.

In the unlikely event that you experience any problems with your unit which cannot easily be remedied, you should contact your supplier and return the unit if necessary. Please include any details of the fault observed and remember to return the unit in its original packing. Techne accept no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier. See the Decontamination Certificate supplied with your unit.

1. Cleaning

Before cleaning your unit ALWAYS disconnect it from the power supply and allow it to cool below 50° C.

Your unit can be cleaned by wiping with a damp soapy cloth. Care should be exercised to prevent water from running inside the unit. Do not use abrasive cleaners.

2. Overtemperature cutout

In the event of no heater power, check the mains plug and lead. Repeated operation of the cutout indicates a serious fault: you may need to return the unit to your supplier for repair.

3. Fuses

Your unit is protected by one or two fuses. These should only be changed by suitably qualified personnel.

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THE TU-20HT

Make sure you have read this manual carefully before you use the thermoregulator.

The Calibration Bath

Techne supply a range of liquid calibration baths from 5 to 12 litres on which you can fit one of the thermoregulator range of units. The bath inner containers are manufactured from stainless steel, for maximum corrosion resistance. The 5 and 7 litre baths are deep drawn with large easy clean corner radii. The 12 litre bath is of welded construction.

Description

The thermoregulator is designed to fit all standard laboratory baths, especially Techne baths. It will heat, circulate and safely control the temperature of the liquid in the bath within precise limits.

The TU-20HT will control the temperature of the liquid in a suitable bath within the range -40°C to 250°C (see the specification for details). However, temperatures from -40°C to "5°C above ambient" require an additional cooling system such as a Techne Fridge unit.

The instrument consists of the following main parts:

- A base moulding in PPS plastic.
- Pump support legs and pump assembly in Victrex Peek.
- A heater assembly in Stainless Steel.
- A base plate made from stainless steel to which are mounted the motor, overtemperature cut-out, mains switch, fuse holder and PCB assembly.
- A cover made from Noryl plastic to which is fitted the digital controls.

The bath temperature is monitored and controlled by a PRT in conjunction with a 3 term controller.

Overtemperature protection is provided by means of an adjustable cutout which also activates an audible alarm. The pump motor and the transformer are also fitted with thermal fuses.

A portable clamp or a bridge piece are available as alternative ways of fixing the thermoregulator to a bath.

Unpacking


When unpacking the unit, check that the following have been removed from the packing: The TU-20HT unit; a Guarantee card; a De-contamination Certificate.

Within the guarantee period, shown on the Guarantee Card, we undertake to supply replacements free of charge for parts which may on examination prove to be defective, provided that the defect is not the result of misuse, accident or negligence. Any instrument requiring service under this guarantee should be sent to the supplier through whom it was purchased, or, in the case of difficulty, it should be carefully packed in its original packing and consigned, carriage paid, to us. Techne takes no responsibility for returned goods damaged in transit.


Returned goods will not be processed without a Returns Authorisation Number. Call + (44) (0)1223 832401 for a number. On all correspondence, please quote the Serial Number in full and/or the Sales Order Number. **Please write the Returns Number on the outside of any packing.**

Set Up


- 1 Ensure that the bath is set up on a flat level surface.
- 3 Fit the unit securely to the bath using the correct bridge piece for the unit/ bath or a portable clamp, see the list of accessories. THE UNIT MUST ALWAYS BE MOUNTED WITH THE BACK AND THE SWITCH END OUTSIDE THE AREA OF THE BATH. This will reduce the infiltration of hot vapours into the cooling system of the thermoregulator. Ensure that at all times the air inlet and outlet remain clear of obstructions. Free circulation of air inside the unit is essential for proper cooling of the electronics and pump motor.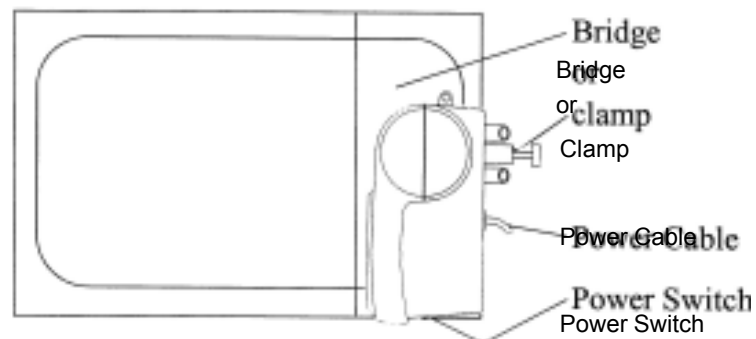


- 3 Use the M4 screw provided with the bridge kit to secure the bridge piece to the thermoregulator at the heater end of the thermoregulator, do not over tighten it. Thumb screw 2 secures both the unit and the bridge to the bath. Thumb screw 3 secures both the cover and the bridge to the bath. Thumb screw 4 is fitted at the Techné and secures the cover to the bridge. A bung is provided to seal off the thermometer hole in the bridge piece if a thermometer is not used. It is necessary to fit this bung to reduce heat losses and prevent steam or other vapours getting to the thermoregulator.

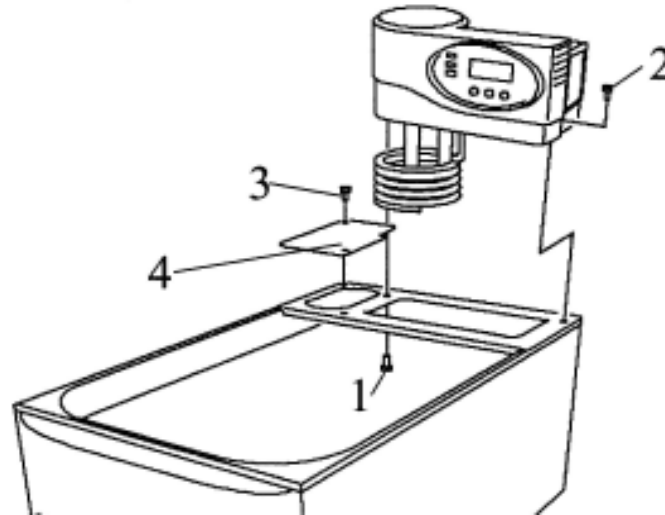- 4 If you do not want to use a bridge piece to secure the unit to a bath then you must use a clamp. The clamp bracket fits between the pump legs in the slot provided. Slide the bracket into the slot and use the two screws in clamp kit to secure the clamp to the unit.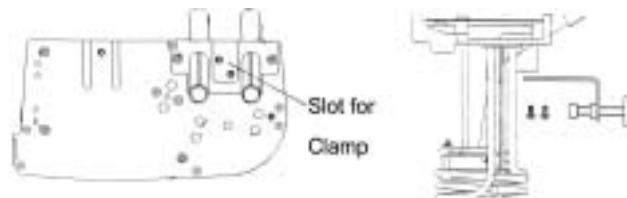

- 5 A cooling coil, which is an optional extra see the list of accessories at the end of this manual, will give control for temperatures between 5°C above ambient and 5°C above the temperature of the water supply.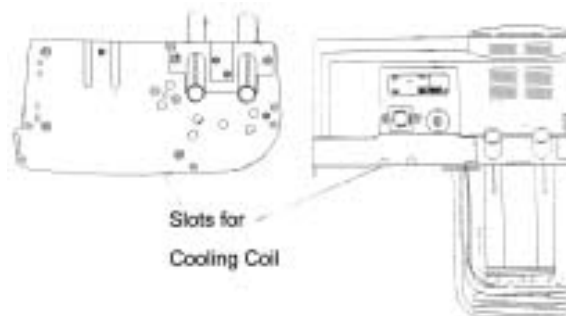


A cooling coil can be fitted to the thermoregulator just below the overtemperature cutout reset button. Connect a hose from a tap to one end of the coil and from the other end of the coil to drain. Adjust the water flow to give the required cooling.

If you are using the thermoregulator on certain baths, for example the Techne Calibration Bath 7 litre, you will not be able to fit a cooling coil. The calibration baths are fitted with their own cooling coil.

- 6 For lower temperatures, a Dip Cooler such as a Techne RU-200 or RU-500 is required. Alternatively you could use either the RB-5A or the RB-12A Techne refrigerated baths, then these have integral fridge coils and no additional cooling is required.

See the instruction Manuals for these units for connections and setup.

- 7 If a thermometer is to be used it may be fitted in the end of the top cover using the O-ring, supplied with the unit, between the two lugs. If a bridge is fitted it is necessary to remove the blanking plug.

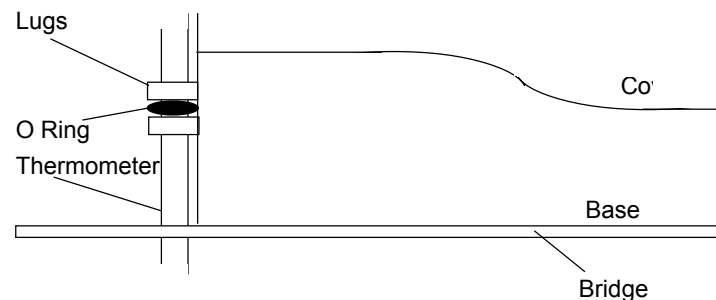- 8 Fill the bath to between the minimum and maximum levels stated in the specification.

If you are using oil at the highest temperatures, remember that they will expand as the temperature rises. Even if you fill the bath with oil to the minimum level, it may well reach the maximum level before the top temperature is reached. Take all precautions and monitor the level through the full range.

If water is used, demineralised water is preferred to reduce the formation of scale. If scale should form, use only mild de-scaling agents to remove it. DO NOT attempt to hammer, chip or scrape the deposits from the surface of the bath

- 9 Recommended liquids:

<i>Temperature</i>	<i>Liquid</i>
-40°C to 0°C	40% water 40% ethylene glycol 20% alcohol
-20°C to 30°C	50% water 50% ethylene glycol
5°C to 95°C	water, preferably de-ionised with neutral pH
10°C to 150°C	Dow Corning Silicone Oil 200 series *
10°C to 250°C	Dow Corning Silicone Oil 210H/100cs series *

* **Warning:** check gel life at top end of range.

Extraction may be necessary at high temperatures; always check the manufacturer's data and safety sheets before using any of the liquids.


In all cases the OVERTEMPERATURE CUT-OUT must be **set correctly** for the liquid being used and the temperature at which it is to operate

- 10 A bath that is fitted with a lid or insulating ball blanket gives the best operating conditions. A lid or ball blanket will prevent vapour loss, heat loss and give better temperature control. If an open bath is used above 80°C (ie where steam or other readily condensing vapours are present) the operation of the unit, particularly the digital display, may be affected.

Below about 80°C a cover becomes less important but will still give better temperature control.

11 The symbols next to the indicator lamps on the front panel of the thermoregulators have the following meanings:

 : the power indicator

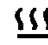 : the heater indicator

 : the overtemperature indicator

12 Symbols on the switch have the following meanings:

I : mains switch On

O : mains switch Off

Specification

The specification was achieved in an 8 litre bath with a ball blanket according to DIN 58966.

Operating temperature range		-40 to +250°C
Working temperature range		5°C above ambient to 250°C
Temperature selection		Digital
Temperature display		Digital LED
Temperature stability		±0.005°C
Set point accuracy		±1°C
Method of control		3 Term
Temperature sensor		PRT
Nominal heater power	230V	1800W
	120V	1500W
	100V	1250W
Maximum watts density		6.2 W/cm ²
Audible alarm for		Overtemperature
Cooling coil		Accessory
Communications		RS232
Protection against hazards		IP30
Safety device classification		2

The TU-20HT has passed:

EN 50081-1:1992

Electromagnetic Compatibility emission standard.

EN 50082-1:1992

Electromagnetic Compatibility immunity standard


(Performance criterion B).

BATHS

Nominal Dimensions

Bath		B-8	B-12	B-18	B-26	B-48
<i>Overall</i>						
Length	A	265 mm	354 mm	530 mm	530 mm	559 mm
Width	B	325 mm	325 mm	325 mm	325 mm	365 mm
Height	C	172 mm	172 mm	172 mm	222 mm	296 mm
<i>Internal</i>						
Maximum Length	D	240 mm	329 mm	505 mm	505 mm	560 mm
Working Length	E	115 mm	205 mm	380 mm	380 mm	435 mm
Width	F	300 mm	300 mm	300 mm	300 mm	330 mm
Max. Working Depth	G	130 mm	130 mm	130 mm	180 mm	255 mm
Min. Working Depth	H	90 mm	90 mm	90 mm	140 mm	215 mm
Maximum Capacity		8 l	12.0 l	18.0 l	26.0 l	48.5 l
Minimum Capacity		6 l	8.4 l	13.2 l	20.5 l	42.5 l

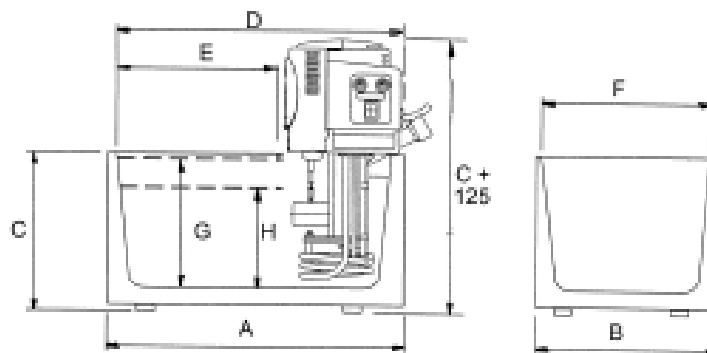**Working Environment**

The TU-20HT units are designed to work safely under the following conditions:

Ambient temp. range 5°C to 40°C

Humidity Up to 95% relative humidity, non-condensing


Note: The control specifications quoted are for an ambient temperature range of 10°C to 30°C. The specification may deteriorate outside this range but the unit will still work safely.

Radio frequency interference tested and passed to EN50081-1.

Immunity Tested and passed to EN50082-1

OPERATION


- Ensure that both the outlet and return pipes have the blanking caps on.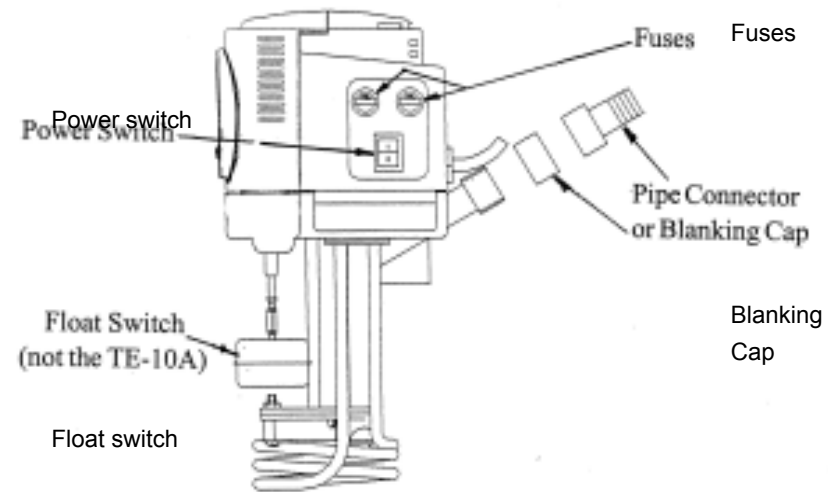

- Switch the unit on by pressing the power switch. The switch and the POWER indicator on the front will light up.

When you Switch On the TU-20HT

When you first switch on, the display will show the edition of the software which your unit uses. For example software issue "5.37" would be shown as follows: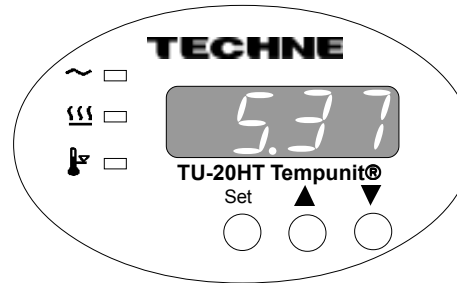


It will display this for 1 second, then the actual temperature of the bath will be indicated.

The Front Panel Controls


The front panel controls consist of three buttons for controlling the display, a four digit LED display and three indicators.

The SET temperature Button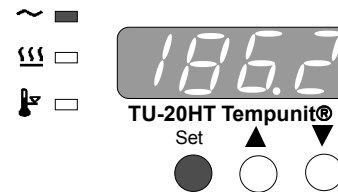


The **SET** temperature button displays the set temperature when pressed.

The UP ARROW Button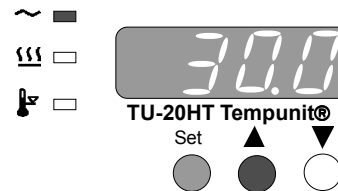

When the **SET** temperature button is held down and the **UP ARROW** button is pressed, the set temperature is increased.

The DOWN ARROW Button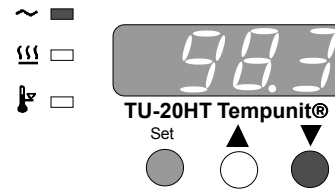


When the **SET** temperature button is held down and the **DOWN ARROW** button is pressed, the set temperature is decreased.

Speed of Change of Set Temperature


Each press of the **UP ARROW** or **DOWN ARROW** buttons will increase or decrease the set temperature by 0.1°C. If the buttons are held down the temperature change will accelerate to 5° per second

Power Indicator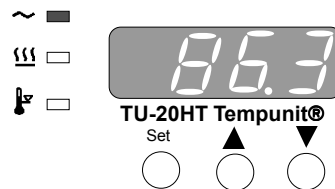


The top indicator shows that there is power to the unit

Heater Indicator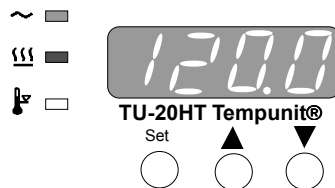


The next indicator shows when the heater is heating. When the temperature is being set, and the new set temperature is higher than the temperature already in the unit, the heater indicator will light as the unit tries to follow the set temperature. If the light is on continuously the heater is getting constant power. The only exception is described under Over-Temperature Indicator. As the temperature approaches the set temperature the heater indicator will flash. When set temperature is reached the indicator will stay on for shorter periods. If the bath temperature is above the set temperature then the indicator will be off, as the heater is not getting any power.

Over-Temperature Indicator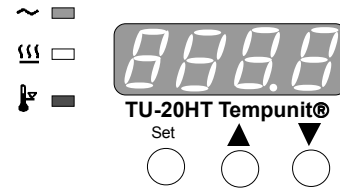


If the unit should, for any reason, exceed the temperature set for the over-temperature cutout, the over-temperature indicator will light. The heater will have been switched off and the unit will begin to cool even if the heater light is on (the light staying on or not depends on which circuit has sensed an over-temperature).

Low Liquid Level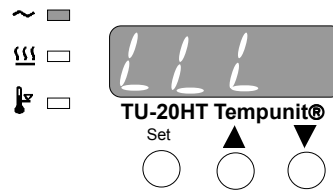

The float switch will trip if the liquid gets below a safe level; the display will change to "LLL". The heater will be switched off. Fill the bath to above the minimum level of 90mm and the display will return to the 'present' temperature; the unit will again work.

Sensor fault Indicator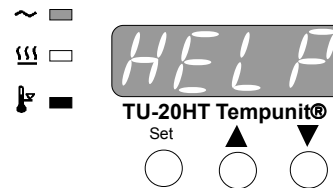

If there should, for any reason, be a sensor fault, the bottom indicator will light. The power to the heater will have been switched off and the unit will begin to return to ambient even if the heater light is on (the light staying on or not depends on which circuit has sensed a fault).

After Use


When you have finished heating samples, remember that parts of the unit and the samples may be very hot. Take the precautions listed earlier. We recommend that the samples should be allowed to cool to 50°C before being removed from the bath. They will still have to be handled with care.

Should you want to remove the unit from the bath, it too should be allowed to cool to 50°C before being removed.

Remember the bridge, the lid (if used), the bath and all other parts close to the bath will be hot while it is in use.

Setting the Overtemperature Cut-out

An adjustable overtemperature cutout is fitted to the TU-20HT. When the cutout operates, the heater will stop working, the audible alarm will sound and the 'Overtemperature cutout indicator' will illuminate. The pump will continue working.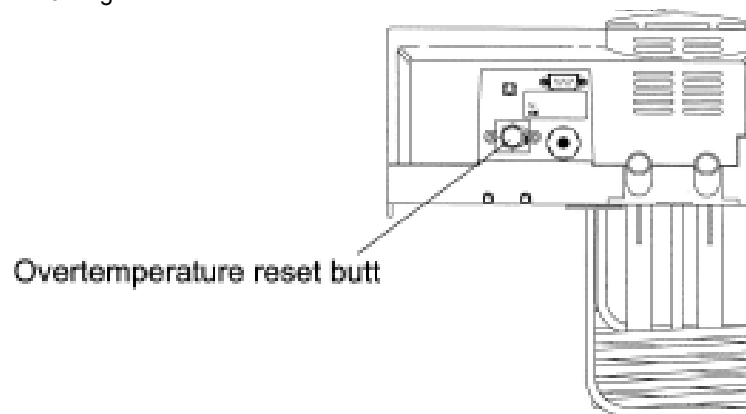

It should be set to approximately 10°C ABOVE THE OPERATING TEMPERATURE. This can be done in one of two ways; '1' is more accurate; '2' is quicker (and better if you do not want to overheat the liquid). For both, first turn the reset button fully clockwise then:

- 1 Heat the bath to the **desired cutout temperature** and turn the reset button anticlockwise until the cutout just trips.

Either: Switch off at the mains power. Remove the unit from the liquid and press the reset button. Return the unit to the liquid, switch on the mains power. The heater will again work.

Or: Allow the liquid to cool, maybe as much as 40°C, and press the reset button. The heater will again work.


- 2 Heat the bath to the required **maximum operating temperature** and turn the reset button anticlockwise until the cutout just trips. Turn the reset button clockwise one small division on the label and press the reset button. The heater will again work.

An alarm will sound when the overtemperature cutout is activated by a temperature higher than that set for the overtemperature cutout. Reset the overtemperature cutout, by pressing the reset button, to stop the alarm.

RS232 SERIAL INTERFACE

The TU-20HT may send data logging information to an IBM PC or compatible computer by connecting the unit and the PC via an RS232 cable, and installing the "Thermsoft" software. Contact your supplier for details.

RS232
Connection

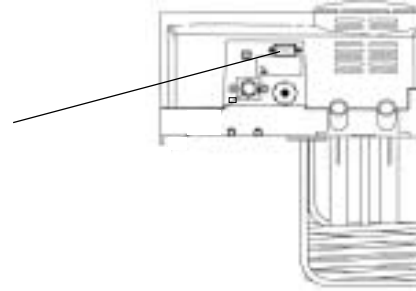The RS232 cable must be fitted to both the unit and the PC before either unit is powered up, otherwise, data integrity cannot be guaranteed. Once the cable is fitted, it does not matter which unit is powered up first.

The following tables indicate the cable specifications for a 9-way PC serial port:

Thermoregulator		PC	
<i>9-way female D type</i>		<i>9-way female D type</i>	
pin	signal	pin	signal
Case	F.GND	Case	F.GND
3	TxD	3	TxD
2	RxD	2	RxD
7	RTS	7	RTS
8	CTS	8	CTS
6	DSR	6	DSR
1	CD	1	CD
4	DTR	4	DTR
5	S.GND	5	S.GND

Sending data to the PC

The procedure for sending data to the PC is described in the "help" instructions as part of the software. The PC must be properly connected by the RS232 cable and running "Thermsoft" to accept or display data.

ADDITIONAL INFORMATION

NOTE THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL. **REMOVING THE TOP CASE EXPOSES POTENTIALLY LETHAL MAINS VOLTAGE.** THERE ARE NO OPERATOR MAINTAINABLE PARTS WITHIN THE EQUIPMENT.

In the unlikely event that you experience any problems with your Thermoregulator which cannot easily be remedied, you should contact your supplier and return the unit if necessary. Please include any details of the fault observed and remember to return the unit in its original packing. Techne accept no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier.

Operator maintenance

1. Cleaning

Before cleaning your unit ALWAYS disconnect from the power supply and allow to cool below 50° C.

Your Thermoregulator can be cleaned by wiping with a damp soapy cloth. Care should be exercised to prevent water from running inside the unit. Do not use abrasive cleaners.

2. Overtemperature cut-out

The overtemperature cut-out is a sensitive mechanical device and mechanical shock can cause it to trip.

- In the event of no heater power, check the mains plug and lead, then reset the cut-out control.
- Repeated operation of the cut-out indicates a serious fault: you may need to return the unit to your supplier for repair.

3. Fuses

Your unit is protected by two fuses.

230V 2 x F10A; 120/100V 2 x T16A

These should only be changed by suitably qualified personnel.

The fuses must only be replaced by a fuse of the same type and value.

If the fuses blow persistently, a serious fault is indicated and you may need to return the unit to your supplier for repair.

Calibration

Remember that if you change the calibration from that set at the factory you may change the calibration at all temperatures. You may get different calibration with different baths and/or liquids.

In order to ensure that the calibration you are setting is correct, you will need to use an independent calibrated probe or thermometer.

Set the temperature display to the particular temperature at which you require to control. Measure the actual temperature of the bath liquid using a calibrated probe or thermometer.

If the calibration is not correct then you can follow this procedure.

- a Hold down the Up and Down buttons and then press the Set button at the same time for 5 seconds.
- b The display will change from the bath temperature to "EEEE".
- c Press the 'SET' button and either the 'UP' or the 'DOWN' button to adjust the display to the same temperature as the measured value.
- d Press 'UP' and 'DOWN' together to confirm the value. The display will return to the bath temperature and the unit will control with the new calibration parameters.

Replacement parts


Each unit is supplied with an O ring for a Thermometer . The following parts may be purchased if replacements or alternatives are required:

<i>Part Number</i>	<i>Description</i>
6007349	O ring (thermometer)
6103913	Cap seal
6103475	Cap

Accessories

There are many accessories both for the thermoregulators and the baths. These are constantly being updated and details can be found in Techne Sales literature. However they do include:

<i>Flat Lids</i>	<i>To fit Bath</i>	<i>Gabled Lids</i>	<i>To fit Bath</i>
FFLAT08	B-8	FGABLE08	B-8
FFLAT12	B-12	FGABLE12	B-12
FFLAT18	B-18 or B-26	FGABLE18	B-18 or B-26
FFLAT48	B-48	FGABLE48	B-48
<i>Bridge</i>		<i>Clamp</i>	
FBRIDGE1	B-8 to B-26 bridge	FCLAMP	TE clamp
FBRIDGE2	B-48 bridge	FCLAMP2	TU clamp
<i>Cooling Coil</i>			
FCC01	Cooling Coil		
<i>Fridge Baths</i>		<i>Dip Coolers</i>	
FRB5D	RB-5A 230V UK	FRU2D	RU-200 230V UK
FRB5E	RB-5A 230V Europe	FRU2E	RU-200 230V Europe
FRB5P	RB-5A 120V	FRU2P	RU-200 120V
FRB2D	RB-12A 230V UK	FRU5D	RU-500 230V UK
FRB2E	RB-12A 230V Europe	FRU5E	RU-500 230V Europe
FRB2P	RB-12A 120V	FRU5P	RU-500 120V
<i>Flow Coolers (not for TU-20HT)</i>			
FFC2D	FC-200 230V UK	Test Tube Racks; and so on.	
FFC2D	FC-200 230V Europe		
FFC2P	FC-200 120V		
FFC5D	FC-500 230V UK		
FFC5D	FC-500 230V Europe		
FFC5P	FC-500 120V		
FTU232	"Thermsoft" kit inc: disc, cable and connector for TU-20HT		

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Thank you for reading this data sheet.

For pricing or for further information, please contact us at our UK Office, using the details below.

UK Office

Keison Products,

P.O. Box 2124, Chelmsford, Essex, CM1 3UP, England.

Tel: +44 (0)1245 600560

Fax: +44 (0)1245 600030

Email: sales@keison.co.uk

Please note - Product designs and specifications are subject to change without notice. The user is responsible for determining the suitability of this product.