
**TC-9D**  
**Temperature**  
**Controller**  
Operator's Manual

Issue 1  
10/20/2011

**CONTENTS****TC-9D Operator's Manual**

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	<i>page</i>
<b>SAFETY and INSTALATION</b>	
English	4
Français	6
Deutsch	8
Español	10
<b>INTRODUCTION</b>	12
Packing	12
Description	12
Specification	13
Installation	14
<b>OPERATION</b>	15
<b>TECHNICAL INFORMATION</b>	16
Operator maintenance	16
Servicing	17
Replacement parts	18
Wiring Diagram	19

**NOTES**

## Introduction

Please read all the information in this booklet before using the unit.

## Warning

HIGH TEMPERATURES ARE DANGEROUS: they can cause serious burns to operators and ignite combustible material.

Techne have taken great care in the design of these units to protect operators from hazards, but Operators should pay attention to the following points:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS;
- DO NOT put hot objects on or near combustible objects;
- DO NOT operate the unit close to inflammable liquids or gases;
- DO NOT place any liquid directly in your unit;
- At all times USE COMMON SENSE.

## Operator Safety

All Operators of Techne equipment must have available the relevant literature needed to ensure their safety.

It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures. If the equipment is used in a manner not specified by Techne the protection provided by the equipment to the Operator may be impaired.

All Techne units have been designed to conform to international safety requirements and are fitted with an overtemperature cutout. On some models, the cutout is adjustable and should be set to suit the application. On all other models the cutout is preset to protect the unit.

If a safety problem should be encountered, switch off at the mains socket and remove the plug from the supply.

## Installation

1. All Techne units are supplied with a power cable. This may be integral or plug-in.
2. Before connecting the mains supply, check the voltage against the rating plate. The rating plate is on the rear of the unit. Connect the mains cable to a suitable plug according to the table below.

**Note that the unit must be earthed to ensure proper electrical safety.**

<i>Connections</i>	<i>220V-240V</i>	<i>110V-120V</i>
Live	Brown	Black
Neutral	Blue	White
Earth	Green/yellow	Green

The fused plug supplied with the mains lead for use in the UK is fitted with the following value fuse to protect the cable: 13AMP.

The fuse in the unit protects the unit and the operator

Note that units marked 230V on the rating plate work at 220V; units marked 120V work at 110V.

In both cases, however, the heating rate will degrade by approximately 8%.

3. Plug the mains cable into the socket on the rear of the unit.
4. Place the unit on a suitable bench or flat workspace, or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction.

## After use

When you have finished heating samples, remember that parts of the unit – the tubes, blocks and associated accessories – may be very hot. Take the precautions listed earlier.

## Guarantee

The unit is guaranteed against any defect in material or workmanship for the period specified on the enclosed guarantee card. This period is from the date of purchase, and within this period all defective parts will be replaced free of charge provided that the defect is not the result of misuse, accident or negligence. Servicing under this guarantee should be obtained from the supplier.

Notwithstanding the description and specification(s) of the units contained in the Operator's Manual, Techne hereby reserves the right to make such changes as it sees fit to the units or to any component of the units.

This Manual has been prepared solely for the convenience of Techne customers and nothing in this Instruction Book shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the units or components.

## Operator maintenance

NOTE: THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL.

REMOVING THE SIDE, FRONT OR REAR PANELS EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES.

THERE ARE NO OPERATOR MAINTAINABLE PARTS WITHIN THE EQUIPMENT.

In the unlikely event that you experience any problems with your unit which cannot easily be remedied, you should contact your supplier and return the unit if necessary. Please include any details of the fault observed and remember to return the unit in its original packing. Techne accept no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier. See the Decontamination Certificate supplied with your unit.

### 1. Cleaning

Before cleaning your unit ALWAYS disconnect it from the power supply and allow it to cool below 50° C.

Your unit can be cleaned by wiping with a damp soapy cloth. Care should be exercised to prevent water from running inside the unit. Do not use abrasive cleaners.

### 2. Fuses

Your unit is protected by one or two fuses. These should only be changed by suitably qualified personnel.

If the fuses blow persistently, a serious fault is indicated and you may need to return the unit to your supplier for repair.

## Contact Information

For technical, sales or servicing information, contact your local Techne dealer or,

Bibby Scientific Ltd  
Beacon Road, Stone  
Staffordshire  
ST15 0SA, United Kingdom  
Telephone: +44(0)1785 812121  
Fax: +44(0)1785 813748  
e-mail: [equipment@barloworld-scientific.com](mailto:equipment@barloworld-scientific.com)

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## INTRODUCTION

Read the whole of this book before commencing work with the unit.

The TC-9D Temperature Controller is designed as an alternative to the energy regulator control provided in the laboratory fluidised baths. It is supplied with a Chromel Alumel thermocouple which fits into the sheath built into the fluidised bath.

The TC-9D provides the following features not provided by the energy regulator:

- 1 Precise temperature setting.
- 2 Improved temperature stability.

Control is achieved by monitoring the signal from the thermocouple sensor. By this means the controller will react directly to correct temperature stability.

- 3 Stable temperature under varying loads.

With the energy regulator, the dial setting must be manually adjusted to correct for changes in heat requirements due to a variation in bath loading or ambient temperature. The TC-9D will automatically adjust the heater output to take account of these varying conditions.

The TC-9D consists of a temperature controller mounted in a case with solid state relay switching and is capable of handling up to 20 amps. Connections to the fluidised bath and to the thermocouple are conveniently made with sockets on the back of the instrument. In use, the TC-9D bypasses completely the energy regulator control on the fluidised bath and cycles all the heating elements on and off at the same time. The unit is mounted on four rubber feet and provided with a three wire cord and thermocouple.

## Packing

When you receive your unit, make sure you keep the original packing in case you ever need to return it for service or repair.

The unit must be transported in the original packing to avoid damage. Techne accepts no responsibility for damage incurred unless the unit is correctly packed and transported in this way.

## Specification

Temperature range	0 to 620°C or 32 to 1148°F
Temperature Stability	±0.2 at 100C (SBL2D) ±0.4 at 100C (SBS4)
Method of control	3 Term PID
Display Resolution	1° digital
Temperature Sensor	Type K
Electrical supply	Any voltage between 85V and 250V.
Overall	
Width	203mm (8")
Depth	305mm (12")
Height	191mm (7.5")
Thermocouple	
Length	152mm (6")
Diameter	5mm (3/16")
Lead Length	2.1m (7ft)

## Installation

- 1 The fluidised bath obtains its power and control through the TC-9D controller. The Fluidised bath has therefore to be connected through the TC-9D to the mains or line supply.

If the fluidised bath has had a plug connected to the lead this needs to be removed. Also remove the Twist lock plug from the rear of the TC-9D. Connect the mains lead from the fluidised bath to this plug.

The connections to the plug are as follows:

Line	Lead	Plug letter
LIVE	Black	Y
NEUTRAL	White	X
GROUND	Green	G

Plug the twist lock plug back into the socket on the TC-9D.

- 2 Plug the thermocouple into the back of the TC-9D.
- 3 Remove the packing bar from the thermostat sheath on the fluidised bath, see the booklet for the bath. It is held in with a nut and olive, a new olive is supplied with the TC-9D. Fit the new olive, as appropriate, to the thermocouple and push the thermocouple fully into the thermostat sheath. Tighten the nut so that the olive holds the thermocouple firmly but do not over tighten.
- 4 Connect a clean dry air supply regulated to 30 PSI **via 3/8"OD** or 9.5mm OD tubing. This tubing should be plugged into the quick disconnect fitting on back of the TC9D labeled "air supply". Connect 3/8" or 9.5mm OD tubing between the Fluidized bath and the quick disconnect fitting on back labeled "Fluidized bath".
- 5 Before connecting the mains supply check the voltage with that marked on the rating plate. Connect the mains cable on the TC-9D to a suitable supply as follows:

CONNECTIONS	230Vor110/120V
Live	Brown or Black
Neutral	Blue or White
Earth	Green/yellow

The rating plate is on the rear of the unit.

- 6 Set the energy regulator on the fluidised bath to 10.
- 7 Switch all the fluidised bath heaters on, as the TC-9D will control the temperature by controlling the power input.

## OPERATION

### TC-9D

- 1 Use the up and down arrows to enter a **setpoint temperature**. The bath will begin to heat immediately. The top display is the measured thermocouple temperature and bottom is the setpoint.
- 2 To change the display from **oC to oF** and vice versa press the 2nd button from the left until the lower display shows "UNITS". Next press Up arrow to reach preferred setting. Press 1st button on the left to exit after changes are made.
- 3 **Two setpoint programming** can be used by entering values at menu option "SP1" and "SP2". See step 2 to access options and exit out. Menu option "SP.SEL" tells the controller which to use.
- 4 A single point **calibration offset** can be applied by accessing option "PV.OFS". See step 2 to access options and exit out. Set a correction value in the opposite direction of your external reference reading. i.e. - reference shows the bath to be 3.0 degrees high, then set "PV.OFS" to "-3.0"

## TECHNICAL INFORMATION

Brief fault finding notes are given in this section.

NOTE THAT ELECTRICAL EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL. REMOVING THE CASE EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES.

THERE ARE NO USER MAINTAINABLE PARTS WITHIN THIS EQUIPMENT.

## User Maintenance

### General

In the unlikely event that you experience any problems with your TC-9D which cannot be easily remedied, you should contact your supplier and return the unit if necessary. Please include details of the fault observed and remember to return the unit in its original packing. Techne accept no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier.

### Fuses

If the temperature controller fails to light when power is supplied to the unit the fuses may have blown. Check that there is no external cause such as a faulty plug.

These should only be changed by a suitably qualified person. If the fuses blow persistently, a serious fault is indicated and you may need to return the unit to your supplier for repair.

### Cleaning your TC-9D

Before cleaning your unit ALWAYS disconnect from the power supply and allow to cool.

You can clean the case of the TC-9D with a cloth dampened in water or ethanol. No part of the unit should be immersed in solvents. Do not use acetone or abrasive cleaners.

## Servicing

Before any servicing of the TC-9D is attempted the unit should be disconnected from the mains.

In the interests of safety both of personal and of the instrument, maintenance and repair work should be carried out only by trained electronic or electro-mechanical technicians. Untrained personnel should not attempt to dismantle the instrument.

### **CONTROLLER**


To remove the controller push the side tabs out to unlock the controller and slide it out of the sleeve. Replacement is the reverse of this. For further details on the controller refer to the Eurotherm operating instructions supplied.

### **SOLID STATE RELAY**

Access is gained to the solid state relay by removing the two screws securing it to the heatsink. Be sure to apply a thin layer of thermal grease when replacing the relay.

### Spare Parts

Item Number	Description	Qty	Notes
1	7002942 COVER	1	
2	7002704 FOOT	4	
3	7002694 TEMPERATURE CONT	1	
4	7002939 TOP LABEL	1	
5	7002941 CHASSIS	1	
6	7002917 THERMOCOUPLE SOCKET	1	
7	7001029 SOCKET MAINS	1	
8	7002915 CABLE CLAMP	1	
9	7002918 MAINS CABLE	2m	
10	7002945 FUSEHOLDER	1	
11	7001395 FUSE 20 amp	1	
12	6100472 HEATSINK	1	
13	7032434 SOLID STATE RELAY	1	
14	7002862 AIR VALVE	1	
15	7002919 FLOW METER (SBS4, SBL1)	1	
16	7002920 FLOW METER (SBL2, SBL2D)	1	
17	7002947 THERMOCOUPLE	1	

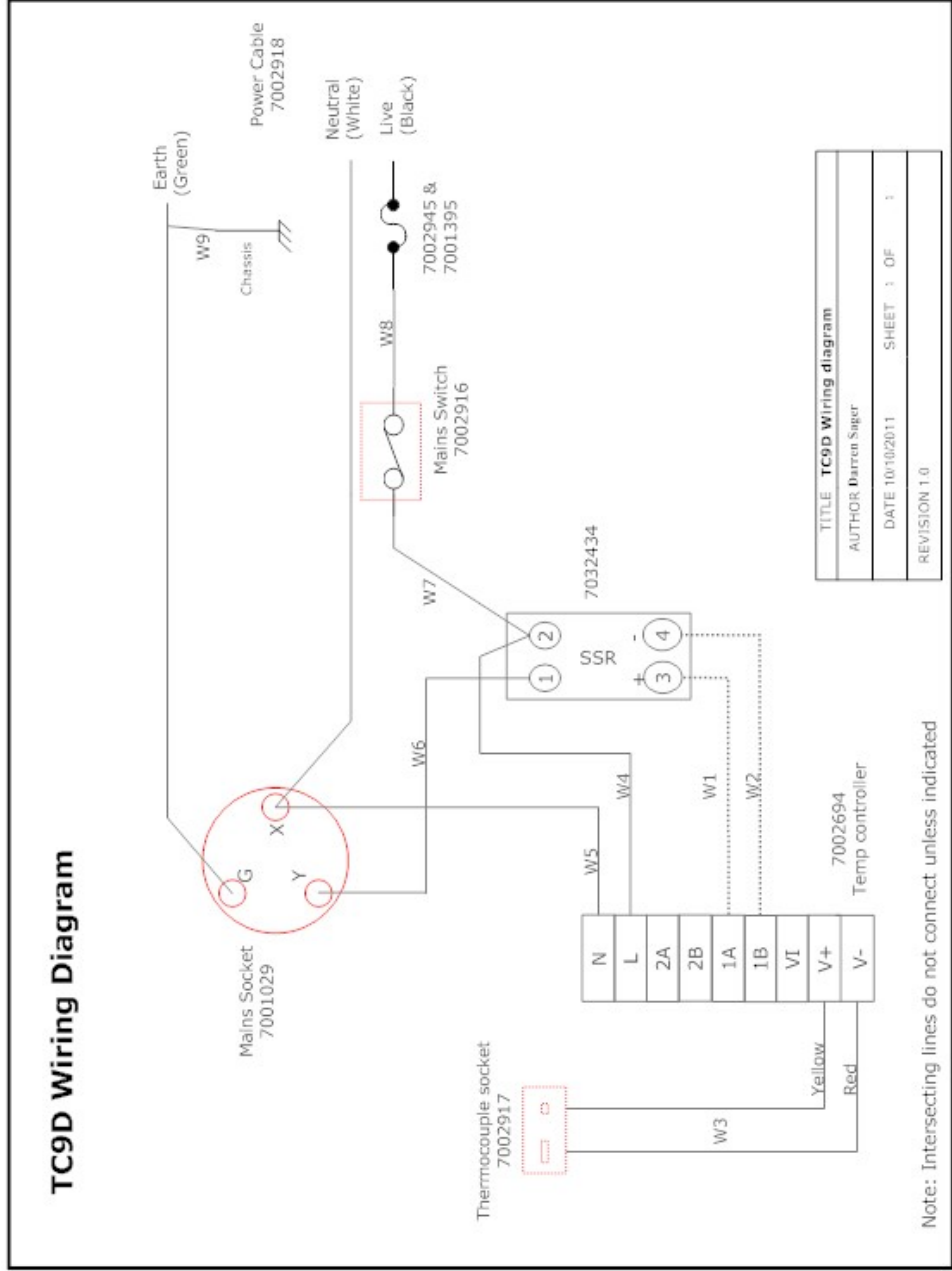Thank you for reading this data sheet.

For pricing or for further information, please contact us at our UK Office, using the details below.

**UK Office**

**Keison Products,**

**P.O. Box 2124, Chelmsford, Essex, CM1 3UP, England.**

**Tel: +44 (0)330 088 0560**

**Fax: +44 (0)1245 808399**

**Email: [sales@keison.co.uk](mailto:sales@keison.co.uk)**

Please note - Product designs and specifications are subject to change without notice. The user is responsible for determining the suitability of this product.